QUICK START GUIDE

ba77037y01 04/2014

HandyLab 100

pH METER

SI Analytics
a xylem brand

Copyright

© 2014, SI Analytics GmbH

Reprinting - even as excerpts - is only allowed with the explicit written authorization of SI Analytics GmbH, Mainz.

Printed in Germany.

HandyLab 100 - Contents

EN English	2
DE Deutsch.....	4
CS Český	6
DA Dansk	8
EL Ελληνικά	10
ES Español.....	12
FI Suomi	14
FR Français	16
HU Magyar	18
IT Italiano	20
JA 日本語	22
KO 한국어.....	24
NL Nederlands	26
NO Norsk	28
PL Polski.....	30
PT Português	32
RO Română	34
RU Русский	36
SK Slovenský	38
SV Svenska	40
TR Türkçe	42
ZH 中文.....	44

HandyLab 100

Overview




The HandyLab 100 meter enables you to perform measurements (pH) quickly and reliably.

The HandyLab 100 meter provides a maximum degree of operating comfort, reliability and measuring certainty.


Keypad

Keys are indicated by brackets <..>.


The key symbol (e.g. <ENTER>) generally indicates a short keystroke (under 2 sec). A long keystroke (approx. 2 sec) is indicated by the underscore behind the key symbol (e.g. <ENTER__>).

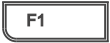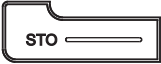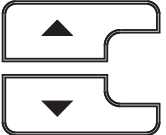	<F1>: <F1__>:	Situation dependent function (is shown on the display) e.g.: <F1>/[Menu]: Opens the menu for measurement settings <F1__>/[Menu]: Opens the menu for system settings
	<On/Off>:	Switches the meter on or off
	<M>:	Selects the measured parameter
	<CAL>: <CAL__>:	Calls up the calibration procedure Displays the calibration data
	<STO>:	Saves a measured value manually
	<RCL>:	Displays the manually stored measured values
	<▲><▼>:	Menu control, navigation
	<ENTER>: <ENTER__>:	Opens the menu for measurement settings / confirms entries Opens the menu for system settings
	<AR>:	Freezes the measured value (HOLD function) Switches the AutoRead measurement on or off

Display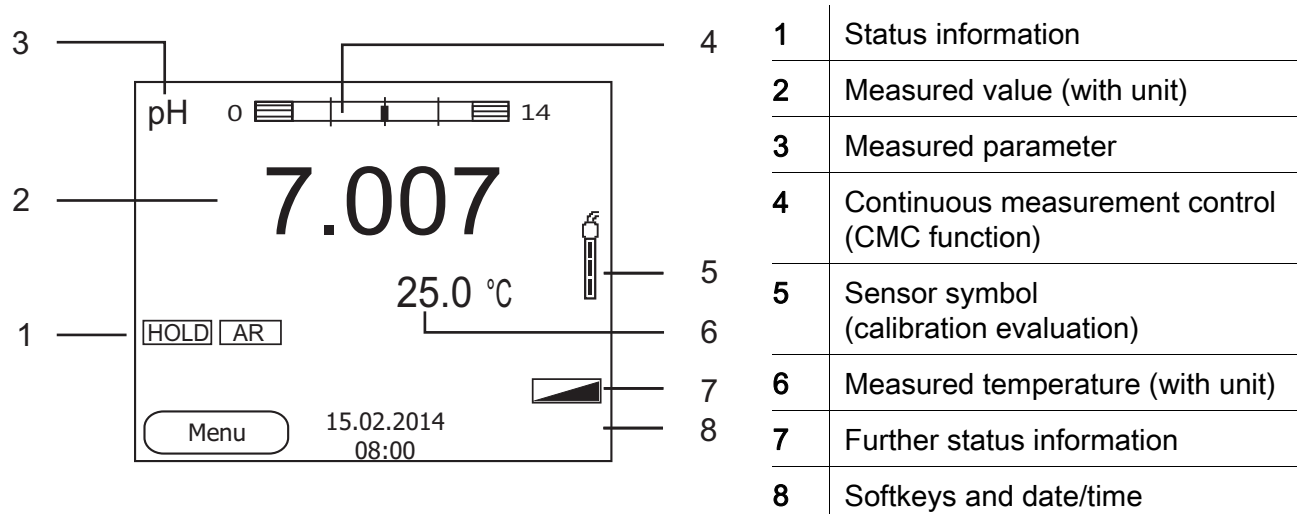

Jack field /connections

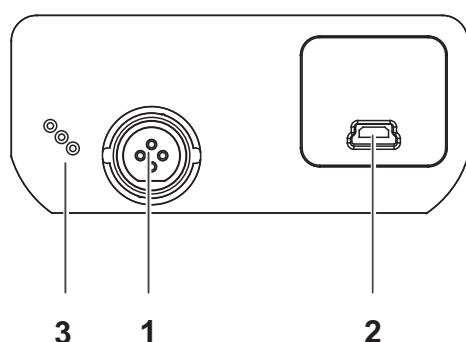1	pH electrode
2	Reference electrode
3	Temperature sensor
4	Service interface

Initial commissioning

- 1 Insert the supplied batteries.
Make sure that the poles are in the correct position.
- 2 Switch on the meter with the <On/Off> key.

Caution

Follow the safety instructions of the used sensors.
The operating manuals of the sensors are available on the supplied CD.

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What can Xylem do for you?

We're 12,500 people unified in a common purpose: creating innovative solutions to meet our world's water needs. Developing new technologies that will improve the way water is used, conserved, and re-used in the future is central to our work. We move, treat, analyze, and return water to the environment, and we help people use water efficiently, in their homes, buildings, factories and farms. In more than 150 countries, we have strong, long-standing relationships with customers who know us for our powerful combination of leading product brands and applications expertise, backed by a legacy of innovation.

For more information on how Xylem can help you, go to www.xyleminc.com

SI Analytics

a **xylem** brand

SI Analytics GmbH

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Tel: +49 (0)6131 / 66 5111

Fax: +49 (0)6131 / 66 5001

D-55122 Mainz

Deutschland, Germany, Allemagne, Alemania

E-Mail: si-analytics@xyleminc.com

www.si-analytics.com

SI Analytics is a trademark of Xylem Inc. or one of its subsidiaries.

© 2015 Xylem, Inc.