QUICK START GUIDE

ba77148y01 10/2015

Lab 855

SI Analytics
a xylem brand

Contents

EN English	3
DE Deutsch.....	7
CS Český	11
DA Dansk	13
EL Ελληνικά	15
ES Español.....	17
FI Suomi	21
FR Français	23
HU Magyar	27
IT Italiano	29
JA 日本語	33
KO 한국어.....	35
NL Nederlands	37
NO Norsk	39
PO Polski	41
PT Português	43
RO Română	47
RU Русский	49
SK Slovensky	51
SV Svenska	53
TR Türkçe	55
ZH 中文.....	57

Copyright

© 2015, SI Analytics GmbH

Reprinting - even as excerpts - is only allowed with the explicit written authorization of SI Analytics GmbH, Mainz.

Printed in Germany.

pH meter, Lab 855

Overview

The compact Lab 855 precision pH meter enables you to perform pH measurements rapidly and reliably.

The Lab 855 handheld meter provides the maximum degree of operating comfort, reliability and measuring certainty for all applications.

The proven calibration procedures and automatic stability control function (AR) support your work with the pH meter.


Keypad

Keys are indicated by brackets <..>.

The key symbol (e.g. <ENTER>) generally indicates a short keystroke (under 2 sec) in this operating manual. A long keystroke (approx. 2 sec) is indicated by the underscore behind the key symbol (e.g. <ENTER_>).


	<On/Off>:	Switches the meter on/off
	<On/Off_>:	Resets calibration data
	<M>:	Selects the measured parameter
	<M_>:	Opens the setting menu for calibration and measurements
	<CAL>:	Calls up the calibration procedure
	<CAL_>:	Displays the calibration data
	<▲>:	Increments values, scrolls
	<▼>:	Decrements values, scrolls
	<ENTER>:	Confirms entries
	<ENTER_>:	Opens the setting menu for system settings

Display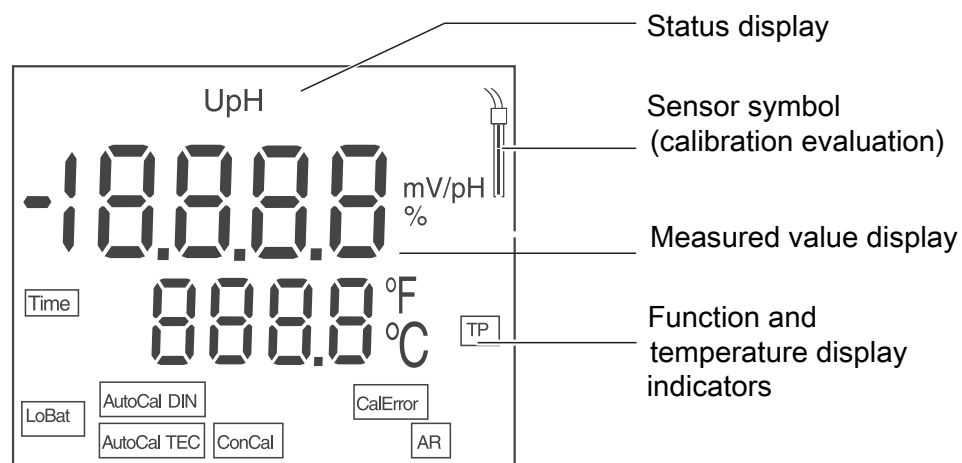

Status display indicators	AutoCal TEC	Calibration with automatic buffer recognition (buffer set: Technical buffers)
	AutoCal DIN	Calibration with automatic buffer recognition (buffer set: DIN buffers)
	ConCal	Calibration with any buffers
	CalError	An error occurred during calibration
	LoBat	With battery operation: Batteries almost empty
	AR	Stability control (AutoRead) is active
	TP	Temperature measurement active
	TIME	Setting of calibration interval

Socket field connections

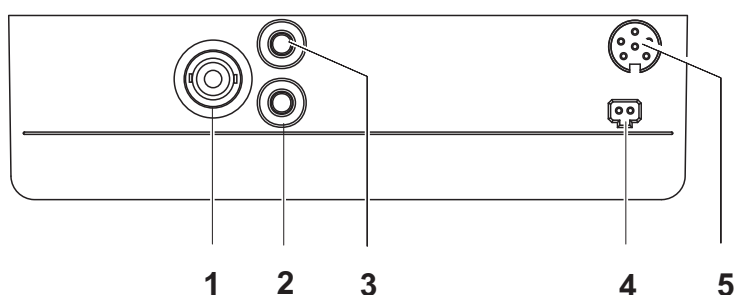1	pH combination electrode
2	Reference electrode
3	Temperature sensor
4	Connection for line adaptor
5	Service interface

Initial commissioning

1	Insert the supplied batteries. When doing so make sure that the poles of the batteries are positioned correctly.
2	Connect the power pack if necessary.
3	Switch on the meter with the <On/Off> key.

Caution

Follow the safety instructions of the used sensors.

The operating manuals of the sensors are available on the supplied CD and on the Internet under www.WTW.com.

Automatic calibration (AutoCal TEC, AutoCal DIN)

1	Press <CAL> repeatedly until the <i>[AutoCal TEC]</i> or <i>[AutoCal DIN]</i> function indicator appears.
2	Confirm the selected calibration procedure with <ENTER> . <i>ct1 [AutoCal TEC]</i> or <i>cd1 [AutoCal DIN]</i> and the temperature are displayed.
3	If necessary, set the temperature of the buffer solution with <▲><▼> .
4	Immerse the pH combination electrode in the first buffer solution.
5	Start the measurement with <ENTER> . The <i>[AR]</i> display indicator flashes. The electrode voltage (mV) or the nominal value of the buffer is displayed.
6	As soon as a stable value is recognized, <i>ct2</i> or <i>cd2</i> appears.

Here you can cancel the calibration procedure with **<M>**. This corresponds to a **single-point calibration**. The value of the zero point (ASY) is displayed. Pressing **<ENTER>** displays the value of the slope (SLO).

7	Thoroughly rinse the combination electrode with distilled water.
8	If necessary, set the temperature of the second buffer solution with <▲><▼> .
9	Immerse the combination electrode in the second buffer solution.
10	Press the <ENTER> key. The <i>[AR]</i> display indicator flashes. The electrode voltage (mV) or the nominal value of the buffer is displayed.
11	As soon as a stable value is recognized, <i>ct3</i> oder <i>cd3</i> appears.

Here you can cancel the calibration procedure with **<M>**. This corresponds to a **two-point calibration**. The value of the zero point (ASY) is displayed. Pressing **<ENTER>** displays the value of the slope (SLO).

12	Thoroughly rinse the combination electrode with distilled water.
13	If necessary, set the temperature of the third buffer solution with <▲><▼> .
14	Immerse the combination electrode in the third buffer solution.
15	Press the <ENTER> key. The <i>[AR]</i> display indicator flashes. The electrode voltage (mV) or the nominal value of the buffer is displayed.

16	The asymmetry is displayed as soon as a stable value is recognized.
17	Press the <ENTER> key. The value of the slope (mV/pH) appears on the display.
18	To return to the measuring mode: Press the <ENTER> key.

Measuring the pH value

1	Immerse the pH combination electrode in the test sample.
2	Using <M> , scroll as necessary until the measured parameter <i>pH</i> is displayed.
3	Wait for a stable measured value. The [AR] display indicator flashes as long as the measured value is not yet stable.

Calibration interval (*Int.C*)

The calibration interval reminds you to calibrate regularly. After the specified calibration interval (*Int.C*) has expired, the symbol for the calibration evaluation flashes. It is still possible to measure.

1	Open the menu for measurement settings with <M__> .
2	Confirm all settings with <ENTER> until <i>Int.C</i> is displayed.
3	Set the calibration interval with <▲><▼> .
4	Confirm the setting with <ENTER> .

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What can Xylem do for you?

We're 12,500 people unified in a common purpose: creating innovative solutions to meet our world's water needs. Developing new technologies that will improve the way water is used, conserved, and re-used in the future is central to our work. We move, treat, analyze, and return water to the environment, and we help people use water efficiently, in their homes, buildings, factories and farms. In more than 150 countries, we have strong, long-standing relationships with customers who know us for our powerful combination of leading product brands and applications expertise, backed by a legacy of innovation.

For more information on how Xylem can help you, go to www.xyleminc.com

SI Analytics

a **xylem** brand

SI Analytics GmbH

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Tel: +49 (0)6131 / 66 5111

Fax: +49 (0)6131 / 66 5001

D-55122 Mainz

Deutschland, Germany, Allemagne, Alemania

E-Mail: si-analytics@xyleminc.com

www.si-analytics.com

SI Analytics is a trademark of Xylem Inc. or one of its subsidiaries.

© 2015 Xylem, Inc.