---

# LF 313T IDS

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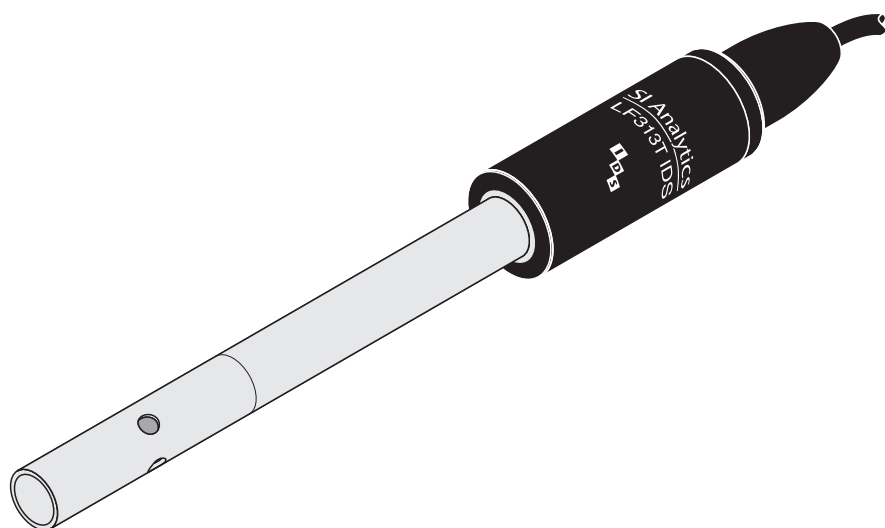
---

# LF 313T IDS

CONDUCTIVITY CELL FOR ULTRAPURE WATER

---

**Copyright**

© 2015, SI Analytics GmbH

Reprinting - even as excerpts - is only allowed with the explicit written authorization of SI Analytics GmbH, Mainz.

Printed in Germany.

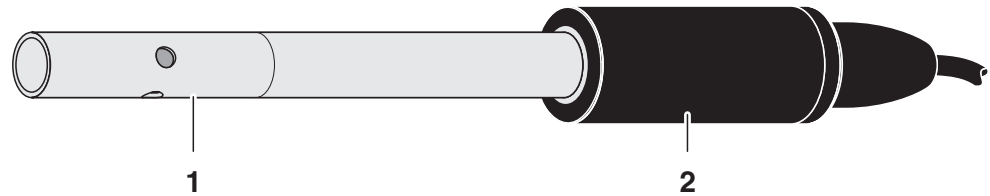
## LF 313T IDS - Contents

<b>1</b>	<b>Overview</b> .....	<b>4</b>
1.1	Structure and function .....	4
1.2	Recommended fields of application .....	4
<b>2</b>	<b>Cleaning</b> .....	<b>5</b>
<b>3</b>	<b>What to do if...</b> .....	<b>5</b>
<b>4</b>	<b>Technical data</b> .....	<b>6</b>
4.1	General data .....	6
4.2	Measuring ranges and resolution .....	7
4.3	Accuracy of the IDS measuring technique .....	8

# 1 Overview

## 1.1 Structure and function

Structure

1	Shaft
2	Connecting head with active electronics

### Automatic sensor recognition

The sensor electronics with the stored sensor data is in the connecting head. The data include, among other things, the sensor type and series number. With each calibration, the calibration data is written in the sensor and the calibration history is recorded. The data is recalled by the meter when the sensor is connected and is used for measurement and for measured value documentation. Storing the calibration data in the sensor ensures that the correct cell constant is automatically used if the sensor is operated with several meters.

The digital transmission technique guarantees the failure-free communication with the meter even with long connection cables. If the sensor firmware is enhanced by SI Analytics, it can be updated via the meter.

## 1.2 Recommended fields of application

Measurements in ultrapure water.

## 2 Cleaning

### Outside cleaning

#### CAUTION

To clean the sensor, disconnect it from the instrument.

We recommend to clean the sensor thoroughly, especially before measuring low conductivity values.

Contamination	Cleaning procedure
Lime sediments	Immerse in acetic acid for 5 minutes (volume share = 10 %)
Fat/oil	Clean with warm water that contains washing-up liquid

After cleaning, thoroughly rinse with deionized water and recalibrate if necessary.

### Aging of the conductivity measuring cell

Normally, the conductivity measuring cell does not age. Special measuring media (e.g. strong acids and bases, organic solvents) or temperatures that are too high may considerably reduce its lifetime or lead to damage. The warranty does not cover failure caused by measuring conditions and mechanical damage.

### Disposal

We recommend to dispose of the measuring cell as electronic waste.

## 3 What to do if...

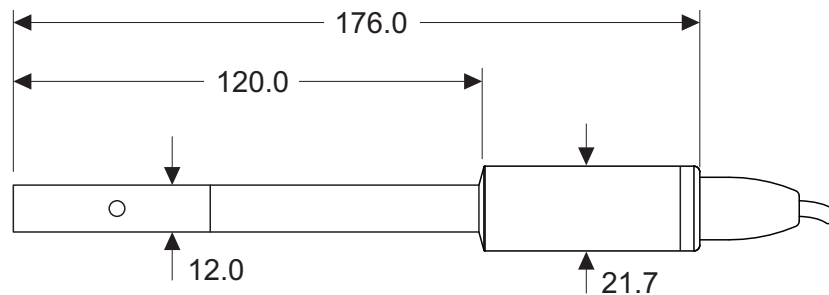
Error symptom	Cause	Remedy
No temperature or conductivity display	<ul style="list-style-type: none"> <li>– No connection between meter and conductivity measuring cell</li> <li>– Cable defective</li> </ul>	<ul style="list-style-type: none"> <li>– Establish connection between meter and conductivity measuring cell</li> </ul>
Measurement delivers implausible conductivity values	<ul style="list-style-type: none"> <li>– Measuring range exceeded</li> <li>– Contamination in the area of the electrodes</li> <li>– Electrodes damaged</li> </ul>	<ul style="list-style-type: none"> <li>– Make sure the correct sensor is being used for the application</li> <li>– Clean the conductivity measuring cell (see section 2).</li> <li>– Return the sensor</li> </ul>
Incorrect temperature display	<ul style="list-style-type: none"> <li>– The temperature sensor is not immersed deep enough in the measuring solution</li> <li>– Temperature sensor defective</li> </ul>	<ul style="list-style-type: none"> <li>– Observe the minimum immersion depth</li> <li>– Return the conductivity measuring cell</li> </ul>

## 4 Technical data

### 4.1 General data

<b>General features</b>	Measuring principle	2-electrodes measurement
	Cell constant	0.100 cm <sup>-1</sup> ± 2 %
	Temperature sensor	Integrated NTC 30 (30 kΩ at 25 °C / 77 °F)

**Dimensions  
(in mm)**

**Weight**      Approx. 90 g (without cable)

<b>Materials</b>	Shaft	Stainless steel 1.4571
	Connection head	POM
	Conductivity electrodes	Stainless steel 1.4571
	Thermistor enclosure	Stainless steel 1.4571

<b>Connection cable</b>	Lengths	1.5 m
	Diameter	4.3 mm
	Smallest allowed bend radius	Fixed installation: 20 mm Flexible use: 60 mm
	Plug type	Socket, 4 pins

<b>Pressure resistance</b>	Sensor with connection cable	IP 68 (2 x 10 <sup>5</sup> Pa or 2 bar)
	Cable plug	IP 67 (when plugged in)

The LF 313T IDS meets the requirements according to article 3(3) of the directive, 97/23/EC ("pressure equipment directive").

<b>Measurement conditions</b>	Conductivity measuring range	0.01 $\mu\text{S/cm}$ ... 200 $\mu\text{S/cm}$
	Temperature range	-5 ... 70 °C (100 °C) 23 ... 158 °F (212 °F)
	Max. admissible overpressure	$2 \times 10^5$ Pa (2 bar)
	Minimum depth of immersion	30 mm
	Maximum depth of immersion (at temperature)	Whole sensor + cable up to 70 °C (158 °F) Sensor shaft only (=120 mm) up to 100 °C (212 °F)
	Operating position	Any
	<b>Storage conditions</b>	Recommended storing method
Storage temperature		0 ... 50 °C (32 ... 122 °F)
<b>Characteristic data on delivery</b>	Temperature responding behavior	$t_{99}$ (99 % of the final value display after) < 20 s
	Accuracy of the temperature sensor	$\pm 0.2$ K

## 4.2 Measuring ranges and resolution

<b>Measured parameter</b>	<b>Measuring range</b>	<b>Resolution</b>
$\kappa$ [ $\mu\text{S/cm}$ ]	0.01 ... 19.99	0.01
	0.0 ... 199.9	0.1
$\rho$ (Resistivity) [ $\text{k}\Omega\cdot\text{cm}$ ]	5.00 ... 19.99	0.01
	20.0 ... 199.9	0.1
	200 ... 1999	1
$\rho$ (Resistivity) [ $\text{M}\Omega\cdot\text{cm}$ ]	2.00 ... 19.99	0.01
	20.0 ... 199.9	0.1
T [°C]	- 5.0 ... + 100.0	0.1

### 4.3 Accuracy of the IDS measuring technique

Measured parameter	Accuracy ( $\pm 1$ digit)
$\alpha, \rho$	$\pm 0.5$ % of measured value
T [°C]	$\pm 0.1$

# What can Xylem do for you?

We're 12,500 people unified in a common purpose: creating innovative solutions to meet our world's water needs. Developing new technologies that will improve the way water is used, conserved, and re-used in the future is central to our work. We move, treat, analyze, and return water to the environment, and we help people use water efficiently, in their homes, buildings, factories and farms. In more than 150 countries, we have strong, long-standing relationships with customers who know us for our powerful combination of leading product brands and applications expertise, backed by a legacy of innovation.

**For more information on how Xylem can help you, go to [www.xylem.com](http://www.xylem.com)**

## SI Analytics

a xylem brand

### **SI Analytics GmbH**

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Tel: +49 (0)6131 / 66 5111

Fax: +49 (0)6131 / 66 5001

D-55122 Mainz

Deutschland, Germany, Allemagne, Alemania

E-Mail: [si-analytics@xylem.com](mailto:si-analytics@xylem.com)

[www.si-analytics.com](http://www.si-analytics.com)

SI Analytics is a trademark of Xylem Inc. or one of its subsidiaries.

© 2015 Xylem, Inc.