

## Technische Information

### Technical Information

## Elektrodengefäß Z 453 Storage vessel Z453



Abbildung Gefäß ohne Elektrode/  
picture without electrode

**SI Analytics GmbH**  
Hattenbergstrasse 10  
55122 Mainz  
Germany  
Telefon +49 (0) 6131/66 5119  
Telefax +49 (0) 6131/66 5001  
e-mail: support@si-analytics.com  
[www.si-analytics.com](http://www.si-analytics.com)

**SI Analytics**  
a xylem brand

### Beschreibung

Das Elektrodengefäß Typ Z 453 dient zur Aufbewahrung und Lagerung von Messketten mit einem Durchmesser von 12 mm. Zum Öffnen und Herausnehmen einer Elektrode wird bei Sicht von oben auf die Elektrode der Bajonett-Verschluß entgegen dem Uhrzeigersinn bis zum Anschlag gedreht. Zur Befüllung wird Wässerungslösung Typ L 911 verwendet. Nicht mit destilliertem Wasser befüllen! Die Füllhöhe beträgt mindestens 2,5 cm, das entspricht ca. 3,2 ml Wässerungslösung. Zum Verschließen wird der Bajonett-Verschluß bei Sicht von oben im Uhrzeigersinn bis zum Anschlag gedreht.

### Lagerung

Elektroden sollten zwischen 0 und 40 °C gelagert werden.

### Technische Daten

Material:	PMMA
Verschluß:	Bajonett
Schlüsselweite:	17 mm

Einbau für Elektroden mit Schaftdurchmesser:	12 mm
---	-------

Bei längeren Elektroden oder solche mit Festkabel, das Elektrodengefäß gegen Umfallen sichern!

### Bestelldaten

285123170	Elektrodengefäß Z 453
285138590	Wässerungslösung L 911, 1000ml Glasflasche
285138560	Wässerungslösung L 9114, 250ml Glasflasche

Technische Änderungen vorbehalten.

### Description

The storage vessel Z 453 is for storage of electrodes with 12 mm diameter. For opening unscrew the bayonet closing by turning it anticlockwise with line of sight on top of the electrode head. The vessel is filled with watering solution L911. Do not fill it with distilled water. The filling level should be at least 2.5 cm, this corresponds to 3.2 ml watering solution. For closure the bayonet closing will be turned clockwise with line of sight on top of the electrode head.

### Storage

Electrodes should be stored at a temperature between 0 and 40 °C.

### Technical details

Material:	PMMA
Closing:	Bayonet
Wrench size:	17 mm

For electrodes with shaft:	12 mm
-------------------------------	-------

For longer electrodes or with fixed cable secure the storage vessel against dropping!

### Order number

285123170	Storage vessel Z 453
285138590	Irrigation fluid L 911, 1000ml glass bottle
285138560	Irrigation fluid L 9114, 250ml glass bottle

Subject to technical amendment.