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## Operating Instructions ..... Page 5 ..... 6

**Important notes:** Before initial operation of the Tempering Jacket for viscometer please read and observe carefully the operating instructions. For safety reasons the Viscometer Tempering Jacket may only be used for the purposes described in these present operating instructions.

Please also observe the operating instructions for the units to be connected.

All specifications in this instruction manual are guidance values which are valid at the time of printing. However, for technical or commercial reasons or in the necessity to comply with the statutory stipulations of various countries, SCHOTT Instruments GmbH may perform additions to the Tempering Jacket for viscometer without changing the described properties.

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## Tempering Jacket for viscometer

**Fig. 1**

The tempering Jacket with holding plate for Cannon – Fenske Routine viscometers is no longer available.

**Fig. 2**

Tempering jacket with holding plate for Ubbelohde Viscometer

- 1 Tempering Jacket
- 2 Holding plate
- 3 Silicone O-ring
- 4 Quick-release closure


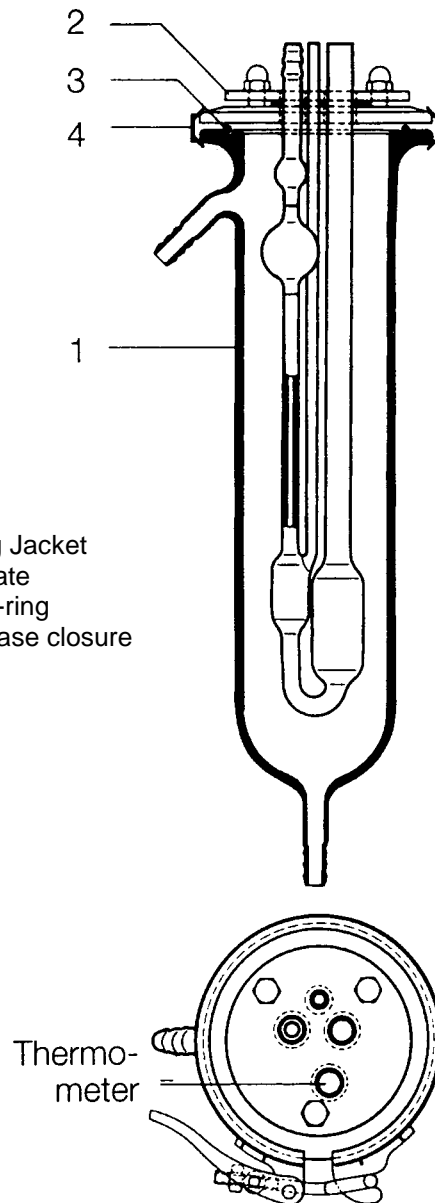
Fig. 3 Explosion drawing Tempering Jacket for viscometer

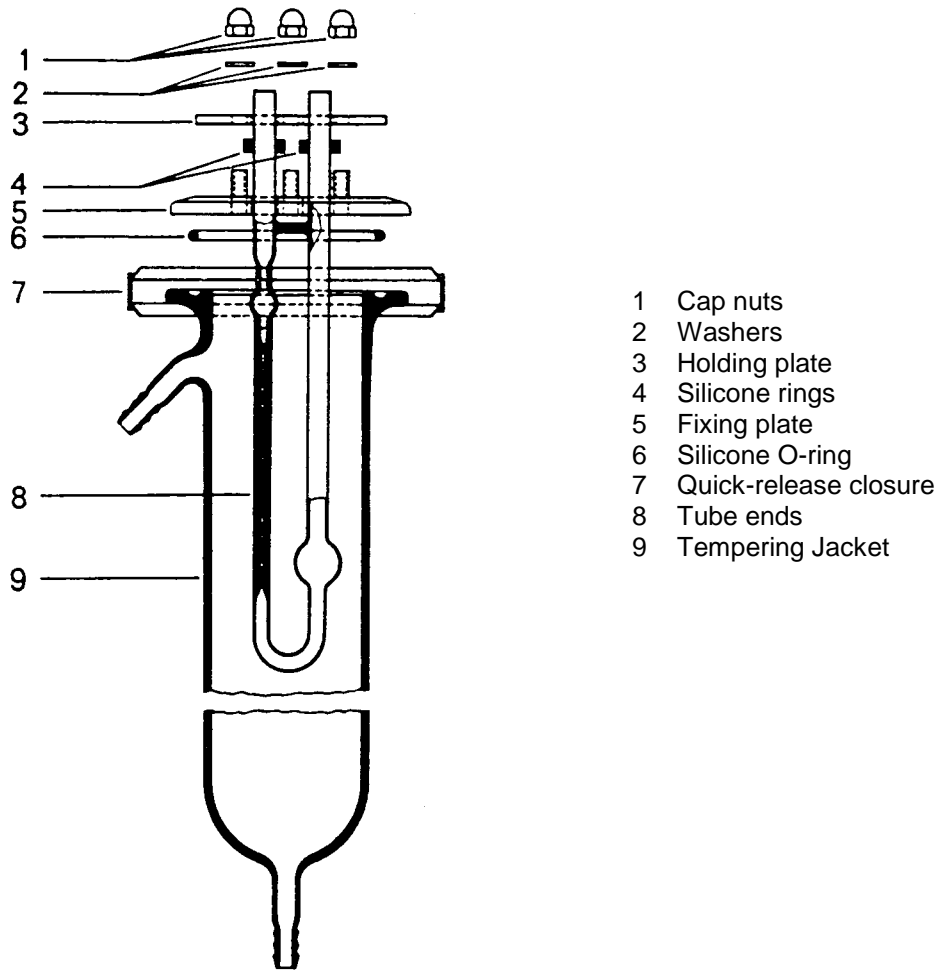Tube ends of viscometer (8) are passed from below through openings in fixing plate (5). Thereafter, silicone rings (4) are slipped over the tube ends and have to be tightened.

The thermometer is inserted in the same way into the aperture provided for this. Holding Plate (3) is subsequently pushed from above over thermometer, tube ends and threaded bolts. By means of cap nuts (1) and washers (2), both plates are pressed against each other under light pressure. Evenly tightening by hand, without using any wrench, normally is sufficient.

O-ring (6) is inserted into groove of the Tempering Jacket (9). Thereafter, the viscometer is inserted into the tempering jacket, and a tight connection is made by means of quick-release closure (7). In order to obtain a tight joint, cap nut on quick-release closure has to be tightened until closure makes contact under light pressure. Level of viscometer is correctly set when upper measuring mark of the viscometer is yet below upper hose connector.

Finally, inlet and outlet are connected to a circulating thermostat by means of flexible tubing (at high temperatures, by means of silicone tubing).

For rinsing, the viscometer is taken out of the Tempering Jacket by opening the quick-release closure.

For measuring and evaluation the current operating instructions for the different viscometers are applied.

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