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**Durchsicht-Thermostat
CT 52**

Operating Instructions


**Transparent Thermostat
CT 52**

Mode d'emploi

**Thermostat transparent
CT 52**

Manual de instrucciones

**Termóstato transparente
CT 52**

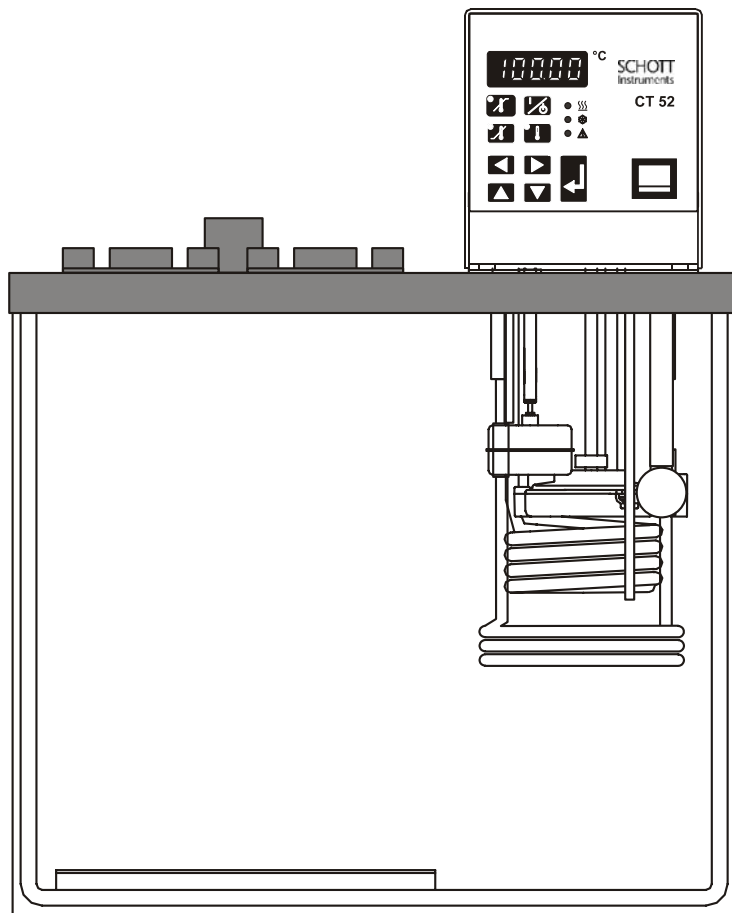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Operating Instructions 27 50

Important notes: Before starting the Transparent Thermostat CT 52 please read the operating instructions carefully. For safety reasons the Transparent Thermostat CT 52 may only be used for the purposes described in this present operating instructions. Please also observe the operating instructions for the units to be connected. All specifications in this instruction manual are guidance values which are valid at the time of printing. However, for technical or commercial reasons or in the necessity to comply with the statutory stipulations of various countries, SCHOTT Instruments GmbH may perform additions to the Transparent Thermostat CT 52 without changing the described properties.

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Supplier's Certificate

We certify that the equipment was verified according DIN EN ISO 9001, part 8.2.4 "Monitoring and measurement of product" and that the specified requirements for the product are met.

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SCHOTT
Instruments

Operating Instructions Transparent Thermostat CT 52

Table of contents

QUALITY ASSURANCE.....	5
YOUR CONTACTS AT SCHOTT INSTRUMENTS GMBH.....	5
SAFETY NOTES	6
DECLARATION OF CONFORMITY	9
1. TECHNICAL DATA OF TRANSPARENT THERMOSTAT CT 52	10
1.1. Technical data immersion circulator CT 52 E	10
1.2. Technical data transparent bath CT 52 B	12
2. OPERATING CONTROLS AND FUNCTIONAL ELEMENTS	13
2.1. Front view CT 52	14
2.2. Rear view CT 52	14
2.3. Overall view	15
3. UNPACKING AND CHECKING	16
4. DESCRIPTION	16
5. PREPARATIONS	17
5.1. Bath liquids	17
5.2. Filling / draining.....	17
6. OPERATING PROCEDURES.....	17
6.1. Power connection	17
6.2. Switching on / Start - Stop	18
6.3. Setting the temperature	18
6.4. Warning functions	19
6.5. Setting the safety temperature (with shutdown function)	20
7. TROUBLESHOOTING GUIDE / ERROR MESSAGES.....	20
8. SAFETY RECOMMENDATIONS	21
9. ATC - ABSOLUTE TEMPERATURE CALIBRATION.....	22
10. ELECTRICAL CONNECTIONS.....	23
11. REMOTE CONTROL	24
11.1. Setup for remote control	24
11.2. Communication with a PC.....	25
11.3. List of commands.....	26
11.4. Status messages	26
11.5. Error messages	27
12. CLEANING THE UNIT	28
13. MAINTENANCE	28
14. DISPOSAL OF THE UNIT.....	28

Quality assurance

Dear customer,

SCHOTT Instruments GmbH implements a **Quality Management System** certified according to EN 29001.

This guarantees the presence of organizational structures which are necessary to ensure that our products are developed, manufactured and managed according to our customers expectations. Internal and external audits are carried out on a regular basis to ensure that our QMS system is fully functional.


Please inform us if, despite our precautionary measures, you should find any product defects. You can thus help us to avoid such faults in future.

Your contacts at SCHOTT Instruments GmbH

Please get in contact with us or the authorized agent who supplied you with the unit if you have any further questions.

SCHOTT Instruments GmbH
Postfach 2443
D-55114 Mainz / Germany
Hattenbergstrasse 10
D-55122 Mainz / Germany

Tel. +49 61 31 / 66 51 11
Fax +49 61 31 / 66 50 01
E-mail: avs@schottinstruments.com
www.schottinstruments.com

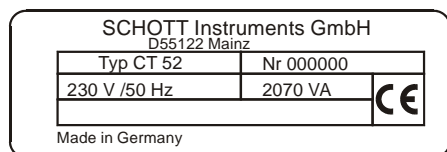The following specifications should be given when product enquiries are made:

- **Unit name** printed on the front of the unit and as specified on the name plate.

Safety notes

These notes are intended to draw your attention to risks which only **you** can recognize and avoid or overcome. They are intended to enhance your own safety consciousness.

We have set the highest quality standards for ourselves and this unit during development and production. Every unit meets relevant safety regulations. **The correct unit usage and proper handling is however solely your responsibility.**

The following notes must be observed:

These operating instructions must be carefully studied! It contains important information on the connection to the local mains supply, correct unit usage and safe handling.

Check for transportation damage during unpacking. Get in contact with supplier and/or carrier for settlement of damage claims. Do not try to start up a damaged unit before the damage has been repaired or you have ascertained the effect of the damage.

Ensure that this manual is always at hand for every unit operator.

Only use this unit solely for the intended application.

Repairs, alterations or modifications must only be carried out by specialist personnel. Considerable damage can be caused by improper repairs. The SCHOTT Instruments GmbH service department is at your disposal for repair work.

Do not operate the unit with wet or oily hands.

Do not expose the unit to spray water or immerse it in water.

Do not clean the unit with solvents (fire risk!) – a wet cloth soaked in household detergent is normally sufficient.

This device is not designed according to the IEC 601-1 standard and should not be operated in rooms used for medical purposes and/or in the vicinity of patients.

**Many unit parts can become hot as a result of normal unit functioning – there is a high risk of burns!
Please ensure that adequate contact protection is provided.**

Do not move the unit from the position where it was set up during operation or when it is still hot. There is a high risk of burns!

Only use the heat transfer liquids recommended by SCHOTT Instruments GmbH . Please refer to the respective EC – Safety Data Sheet.

The temperature controlling i.e. immersing of test tubes, Erlenmeyer flasks or similar objects directly within the circulator constitutes normal circulator practise.

We do not know which substances are contained within these vessels. Many substances are:

- inflammable, easily ignited or explosive
- hazardous to health
- environmentally unsafe

i.e.:dangerous

You alone are responsible for the handling of these substances!

Our advice:

- If in doubt, consult a safety specialist.
- Read the product manufacturer's or supplier's "EC – SAFETY DATA SHEET"
- Read relevant regulations concerning dangerous materials.
- Observe relevant guidelines for laboratories in your country.

The following measures were taken for the protection of the operator:

- Protection Class I according to VDE 0106 T1 (IEC536) i.e. protection against electric shocks by grounding all parts which carry the risk of electric contact.

- This device must only be connected to mains respectacles with a protective ground.

- Protection IP 20 according to EN 60529, i.e. regarding the protection against accidentally touching live parts and damage by foreign matter, it has been ensured that foreign bodies with a thickness or diameter of more than 12 mm cannot penetrate.

No special precautions were taken against the penetration of water and dust. The device should therefore not be used in a dusty atmosphere or in the neighborhood of spray water.

Do not insert wires or tools in any of the openings.

Complete separation from mains is required when:

- all dangers caused by this device are to be avoided,
- cleaning is carried out;
- repairs or maintenance work is about to be carried out

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We declare under our sole
responsibility that the
product

**Transparent
Thermostat
CT 52**

to which this declaration
relates is in conformity with
the standards

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Translation of the legally binding German version

15. Technical data of Transparent Thermostat CT 52

(Version 1 July 2001)

15.1. Technical data immersion circulator CT 52 E

CE sign:

according to the Council Directive 89/336/EEG (electromagnetic compatibility)
generic emission according to standard EN 50 081, part 1
generic immunity according to standard EN 50 082, part 1
according to the Council Directive 73/23/EEG of the (low voltage directive)
last modified by the Council Directive 93/68/EEG
testing basis EN 61 010 part 1.

Country of origin:

Germany / Made in Germany

Temperature ranges:

Operating temperature range with additional cooling [°C]: +10 ... +60
Working temperature range according DIN 58 966
(20 °C ambient temperature) [°C]: +30 ... +60
with tap water cooling* [°C]: +20 ... +60
with flow-through-cooler CK 300 [°C]: +10 ... +60
Temperature accuracy [K]: ±0.02
Temperature drift at changing ambient temperature
between 5 and 40 °C: max.0.002 K/1 °C

Heater capacity:

with 230 V, [VA]: 2000; on request 115 V, [VA]: 1000

Pump:

pressure / flow rate max. [hPa / l/min]: 250/10

Controller:

PID control

Connections:

mechanical:

1. nozzles 2 x 12 mm dia. for flow-through-cooler (CK 300)
2. nozzles 2 x 8 mm dia. for tap water cooling

electrical:

1. connector: control cable to the flow-through-cooler (CK 300)
2. connector: RS232C interface

Data transfer:

Interface: bi-directional serial acc. to EIA RS232C

Data format: selectable

Baudrate: 1200, 2400, 4800 or 9600

Parity: no / odd / even

Length: 7 data-bits

Stop bit: 1

Protocol: CTS/RTS: with or without

Mains power connection:

230 V~, 50 Hz; on request 115 V~, 60 Hz

Technical data of Transparent Thermostat CT 52

Total power consumption: at 230 V, [VA]: 2070; at 115 V, [VA]: 1070

<u>Measurements:</u>	Total, incl. heater, cooling coil and line cord connection: (W x D x H) [cm]: 38 x 26 x 34 Housing above the cover plate of the baths: (W x D x H) [cm]: 12 x 15.5 x 14.5
<u>Weight:</u>	Approx. [kg]: 5
<u>Environment:</u>	Ambient temperature: +10 ... +40 °C (for storage and transportation) Air humidity acc. DIN EN 61 010, part 1: Max. rel. humidity 80 % for temperatures up to +31 °C, linear decrease down to 50 % rel. humidity at a temperature of +40 °C
<u>Safety measures:</u>	Variable overtemperature protection, adjustable from [°C] :+20 .. +80 Low liquid level protection: fixed according to DIN 12 876 safety class III
<u>Protection class:</u>	IP 20 acc. EN 60 529
<u>Power supply:</u>	acc. to class 1, VDE 0106 T1 not for use in explosive atmosphere

* If tap water is used, an operating temperature of appr. 3 °C above the given tap water temperature can be maintained.

15.2. Technical data transparent bath CT 52 B



<u>Filling volume:</u>	[L]: 18
<u>Measurements:</u> W x D x H	[cm]: 38 x 24 x 32
<u>Weight:</u>	[kg]: ca. 4.6, without bath liquid
<u>Bath liquid:</u>	Water
<u>Temperature range:</u>	[°C]: +10 ... +60
<u>Material:</u>	Acryl, high-grade steel
<u>Country of origin:</u>	Germany / Made in Germany

16. Operating controls and functional elements


Front view

- | | | |
|---|---|---------------------------------------|
| 1 |  | Mains power switch, illuminated |
| 2 |  | Start / stop key |
| 3 |  | Working temperature |
| 4 |  | High temperature warning limit |
| 5 |  | Low temperature warning limit |
| 6 |  | MULTI-DISPLAY temperature indication |
| |  | Indicator light - Heating |
| |  | Indicator light - Cooling |
| |  | Indicator light - Alarm |
| 7 |  | Cursors left/right |
| 8 |  | Edit keys (increase/decrease setting) |
| 9 |  | Enter key (store) |


Rear view

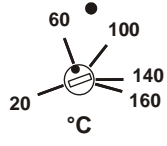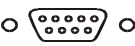- | | | |
|----|---|---|
| 10 |  | Adjustable excess temperature protection
(DIN 12876, safety temperature) |
| 11 | | Threaded fitting (10 mm) for stand rod attachment |
| 12 | | Mains power cable with plug |
| 13 |  | RS232C interface |
| 14 |  | 5-pin connecting socket for cooling unit CK 300 via adapter cable. |


16.1. Front view CT 52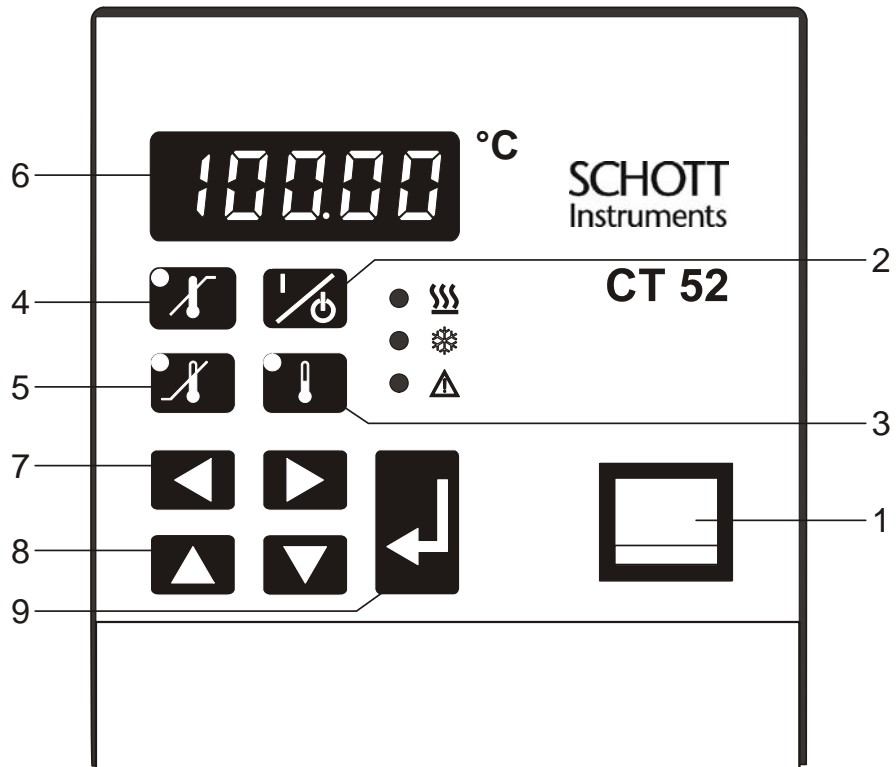

16.2. Rear view CT 52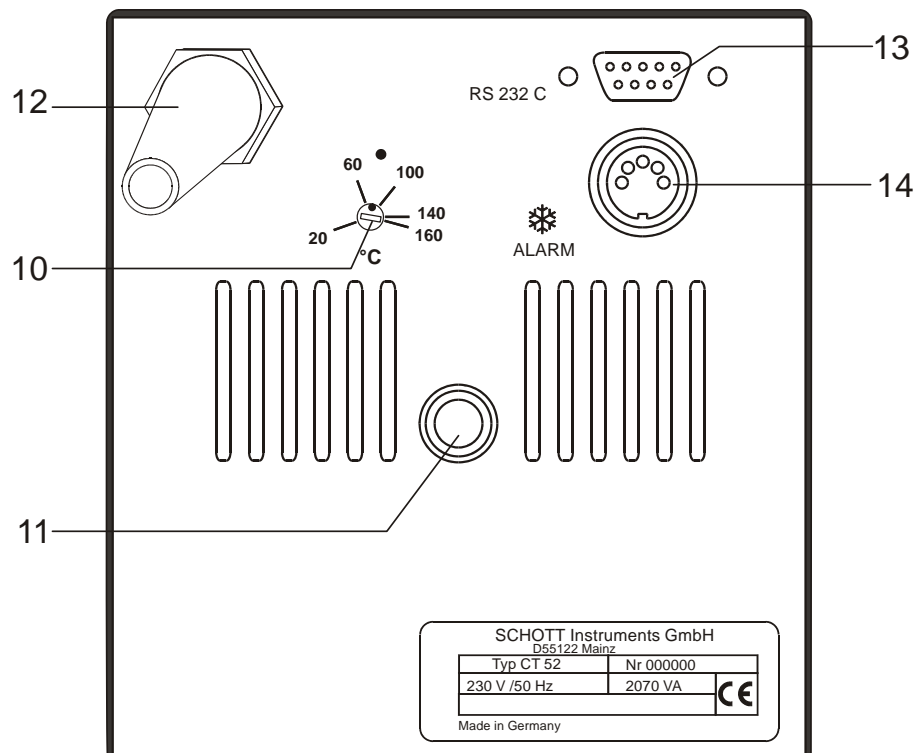

16.3. Overall view

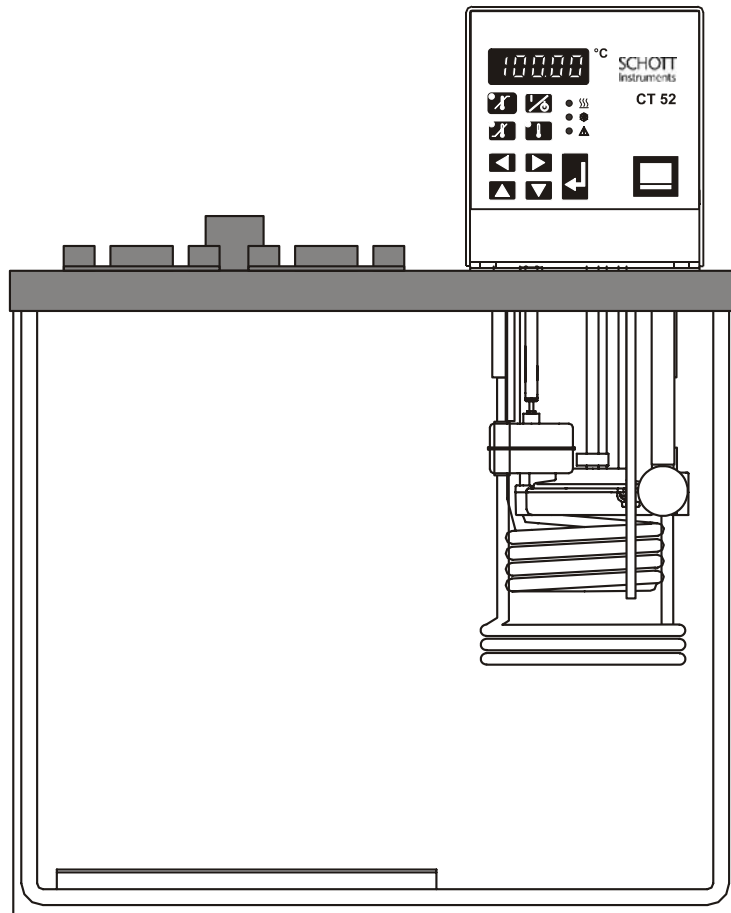Basic unit und transparent bath are supplied ready for operation.

17. Unpacking and checking

Unpack the circulator and accessories and check for damages incurred during transit. These should be reported to the responsible carrier, railway, or postal authority, and a request for a damage report should be made. These instructions must be followed fully for us to guarantee our full support of your claim for protecting against loss from concealed damage. The form required for filing such a claim will be provided by the carrier.

18. Description

The transparent thermostat CT 52 is designed for the heating of water.

The main functional elements are the heater, circulation pump, and control electronics. An electronic proportional temperature control (PID characteristic) adapts the heat supplied to the thermal requirements of the bath. Setting is rapid and simple using the keypad with its easy to learn symbols. Keypad is splash-proof, easily cleaned and ergonomically designed.

The built-in microprocessor enables adjustment and setting of the working temperature and the desired overtemperature and low temperature warning limits. The adjusted settings are displayed on MULTI-DISPLAY.

The safety value for excess temperature protection, a safety installation independent from the control circuit, is adjustable on the rear.

The transparent thermostat CT 52 conforms to the safety requirements specified by DIN 12 876 (safety class III), as well as DIN 58 966, the guideline for first voltage range EN 61010-1 and the NAMUR recommendation.

19. Preparations

19.1. Bath liquids

The transparent thermostat CT 52 is designed for operation with water only!

As a precaution against corrosion, SCHOTT Instruments GmbH recommends the use of Di-water (deionized water) and tap water at a mixture of 1 : 1 to 5 : 1, depending on the hardness of the water.

19.2. Filling / draining

Take care that no liquid enters the interior of the immersion thermostat CT 52 E.

Filling

- Recommended maximum filling level with water as bath liquid is 33 mm below the cover plate.

Draining

- Press the mains switch to turn the transparent thermostat CT 52 off.
- Disconnect the power plug from the mains socket!
- Empty the bath tank (with water dispense pump, for example).

Exercise CAUTION when emptying hot bath liquids!

20. Operating procedures

20.1. Power connection

**Connect the unit only to a grounded mains power socket!
We disclaim all liability for damage caused by incorrect line voltages!**

Check to make sure that the line voltage matches the supply voltage specified on the identification plate.
Deviations of $\pm 10\%$ are permissible.

20.2. Switching on / Start - Stop

Switching on:

Turn on the mains power switch.

The unit performs a self-test. All segments of the 4-digit MULTI-DISPLAY and all indicator lights will illuminate.

Then the software version (example: n 3.03) appears. The display "OFF" or "r OFF" indicates the unit is ready to operate (standby mode).

The circulator enters the operating mode activated before switching the circulator off:

keypad control mode (manual operation) or
remote control mode (operation via personal computer).

Start:

- Press the start/stop key.
- - The MULTI-DISPLAY indicates the actual bath temperature. (example: 21.33 °C)

Stop:

- Press the start/stop key.
The MULTI-DISPLAY indicates the message "OFF".

The unit also enters the safe operating state "OFF" or "r OFF" after a mains power interruption. The temperature values entered via the keypad remain in memory. With the circulator in keypad control mode, press the start/stop key to restart operation. With the circulator in remote control mode, the personal computer must first resend the parameters set via the interface before the circulator may be restarted.

20.3. Setting the temperature

Take care of the limited temperature range for use with plastic bath tanks.
The working temperature range is stated in the technical specifications (see page 10)

Setting the working temperature

- ① Press the setpoint key

The indicator light **flashes** and the value previously set appears on the MULTI-DISPLAY.

- ② Use the cursor keys to move left or right on the display until the numeral you wish to change is blinking.


- ③ Use the edit keys to increase or decrease the numeral value (-, 0, 1, 2, 3, ... 9).

- ④ Press enter to store the value (example: 37.00 °C).

The working temperature is maintained constant after a short heat-up time (e. g. 37.00 °C).

20.4. Warning functions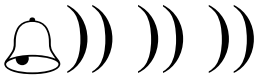

More protection for your samples in the bath!
An audible signal sounds in intervals when the actual temperature value exceeds one of the set limits (patented).

Setting the high temperature limit

- ① Press the key .

The indicator light **flashes** and the value previously set appears on the MULTI-DISPLAY.

- ② Use the cursor keys   to move left or right on the MULTI-DISPLAY until the numeral you wish to change is blinking.

- ③ Use the edit keys   to increase or decrease the numeral value (-, 0, 1, 2, 3, ... 9).

- ④ Press enter  to store the value (example: 39.00 °C).

Setting the low temperature limit:

- ① Press the key .

- ② Follow the instructions


- ③ for .

- ④ (example: 35.00 °C).


Recommendation:

Set the high temperature limit for all applications. The low temperature limit is particularly important for continuous operation.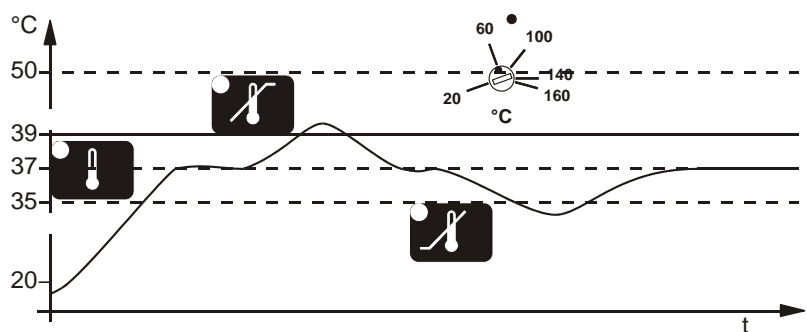

20.5. Setting the safety temperature (with shutdown function)

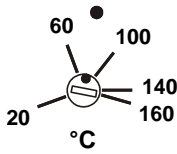(excess temperature protection)

Use a screwdriver to adjust the overtemperature protection device at the rear of the thermostat to the desired value (e.g. 80 °C).

Setting range: 20 °C to 160 °C


This safety installation is independent of the control circuit. When the temperature of the bath liquid has reached the safety temperature, a complete shutdown of the heater and pump is effected.

The alarm is indicated by optical and audible signals (continuous tone) and on the MULTI-DISPLAY appears the error message "Error 01".

Recommendation:

Set the safety temperature at 5 to 10 K above the working temperature setpoint.

21. Troubleshooting guide / Error messages

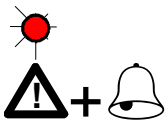Whenever the microprocessor electronics registers a failure, a complete shutdown of the heater and circulation pump is performed.

The alarm light "▲" illuminates and a continuous signal tone sounds.

Cause	Remedy
The circulator is operated without bath liquid, or the liquid level is insufficient.	Replenish the bath tank with the bath liquid.
Tube breakage has occurred (insufficient filling level due to excessive bath liquid pumped out).	Replace the tubing and replenish the bath tank with the bath liquid.
The float is defect (e. g., because damaged in transit).	Repair by SCHOTT Instruments GmbH Hattenbergstrasse 10 D55122 Mainz
The safety temperature value lies below the working temperature setpoint.	Set the safety temperature to a higher value.
A heat reaction or sudden temperature increase, e.g. caused by an exothermic chemical reaction or by the immersion of preheated samples.	Set the safety temperature to a higher value.

Cause

The wires of the working temperature sensor are interrupted or short-circuited.


Remedy

Repair by
SCHOTT Instruments GmbH
Hattenbergstrasse 10
D55122 Mainz

}
|
}
|
}
|
}

other errors

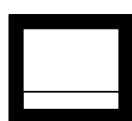After eliminating the malfunction, press the mains power switch off and on again to cancel the alarm state.
If the unit cannot be returned to operation, contact an authorized service station.

Disturbances that are not indicated.

Pump motor overload protection
The pump motor is protected against overloading. After a short cooling interval, the motor will automatically start running.

22. Safety recommendations

Follow the safety recommendations to prevent damage to persons or property. Further, the valid safety instructions for working places must be followed.

- Connect the unit only to a grounded mains power socket
- Ensure secure attachment of heating immersion circulators
- Some parts of the bath cover and the pump connections may become extremely warm during continuous operation. Therefore, exercise particular caution when touching these parts.
- Exercise caution when emptying hot bath liquids!
- Employ suitable connecting tubing.
- Make sure that the tubes are securely attached.
- Avoid sharp bends in the tubing, and maintain a sufficient distance from surrounding walls.
- Regularly check the tubing for material defects (e.g. for cracks).
- Before cleaning the unit, disconnect the power plug from the mains socket.

23. ATC - Absolute Temperature Calibration

37.00

Thermostat (T_T)

36.80

Measuring point (T_M)

Atc0

Atc1

-0.20

37.00

Measuring point T_M

ATC serves to compensate a temperature difference that might occur between circulator and a defined measuring point in the bath tank because of physical properties.

The difference temperature is determined ($\Delta T = T_M - T_T$) and stored as correcting factor (example $\Delta T = -0.2$ K).

- Press the cursor key and enter **at the same time.**
- The MULTI-DISPLAY indicates "Atc0".
- With the edit keys select "Atc1" and then press enter .
- Using the cursor keys and the edit keys set the correcting factor and then press enter (example -0.20 K)
- Press and **at the same time.**

The temperature on the measuring point rises to a temperature of 37.00 °C and is indicated on the MULTI-DISPLAY.

Note:

The correcting always affects the actual working temperature, even if this is set via the interface


The ATC function stays activated until resetting to 00.00 °C.

24. Electrical connections

Connection of circulator CK 300.

The circulator CK 300 is connected to pins 1 and 3. For the connection use the shielded control cable supplied with the circulator.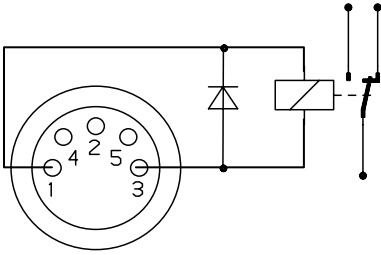

Circuit: Operation = relay powered
 Alarm = relay not powered

Pin assignment:

Pin 1: +24 V (max. current 25 mA)
 Pin 2: 0 V
 Pin 3: Alarm relay
 Pin 4: Reserved - do not use!
 Pin 5: Cooling pulse

RS232C serial interface

This port can be used to connect a computer with an RS232C cable for remote control of the circulator.

Pin assignments:

Pin 2	RxD	Receive Data
Pin 3	TxD	Transmit Data
Pin 5	0 VD	Signal GND
Pin 6	DTR	Data terminal ready
Pin 7	RTS	Request to send
Pin 8	CTS	Clear to send

Interface correspondence:


Circulator (9-pole plug)		Computer (25-pole connector)
Pin 2 RxD	↔	Pin 2 TxD
Pin 3 TxD	↔	Pin 3 RxD
Pin 5 GND	↔	Pin 7 GND
Pin 6 DTR	↔	Pin 6 DSR
Pin 7 RTS	↔	Pin 5 CTS
Pin 8 CTS	↔	Pin 4 RTS

Use shielded cables only.

25. Remote control

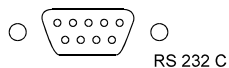25.1. Setup for remote control

RS 232 C

Interface parameters for the circulator are adjusted at configuration level.

Enter or exit the configuration level by pressing the right arrow  and enter  at the same time.

The menu item "Atc" appears on the MULTI-DISPLAY. After pressing the cursor key , the second menu item "REMOTE" is indicated. The interface parameters cannot be adjusted unless the display reads "r 0".

Adjusting interface parameters:

Example: changing the baud rate.

① Using the arrow keys   move the cursor to the desired menu. The actual parameter is displayed (example: "br 24" = 2400 bauds).

② With the up and down arrows   change the number to the desired parameter (example: 4800 bauds). The changed numerals are blinking.

③ The new parameter is stored in memory by pressing enter .

Adjustable interface parameters

REMOTE 0 = keypad control mode
 1 = remote control mode via RS232C

BAUDRATE 12 = 1200 bauds
 24 = 2400 bauds
 48 = 4800 bauds
 96 = 9600 bauds

PARITY 0 = no parity
 1 = odd parity
 2 = even parity

HANDSHAKE 0 = Xon/Xoff protocol (software handshake)
 1 = protocol RTS/CTS (hardware handshake)

Factory settings of the serial interface RS232C:

Baudrate:	4800 bauds
Parity:	2 (even)
Handshake:	1 (hardware handshake)
Data bits:	7
Stop bits:	1

Like all parameters which can be entered through the keypad, interface parameters are stored in memory even after the circulator is turned off.


25.2. Communication with a PC

Suitable terminal programs for communicating with a PC are:

- MS-Windows - TERMINAL.EXE (included with MS-Windows).
- MS-DOS - Procomm Plus, Datastrom Technologies.
- MS-DOS - Norton Utilities.

If the circulator is put into remote control mode via the configuration level, the display will read "r OFF" = REMOTE STOP.

The circulator is now operated via the computer.

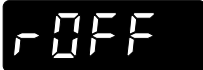In general, the computer (master) sends commands to the circulator (slave). The circulator sends data (including error messages) only when the computer asks for it..

A transfer sequence consists of:

- command
- space (↔; Hex: 20)
- parameter (the character separating decimals in a group is the period)
- end of file (↵; Hex: 0D)

The commands are divided into **in** or **out** commands.

in commands: asking for parameters to be displayed

out commands: setting parameters

The out commands are valid only in remote control mode

Examples:

- Command to set the working temperature to 55.5 °C:
out_sp_00↔55.5↵
- Command to ask for the working temperature :
in_sp_00↵
- Response from the circulator:
55.5↵

25.3. List of commands

Command	Parameter	Response of circulator
version	none	Number of software version (V X.xx)
status	none	Status message, error message (see below)
out_mode_05	0	Stop the circulator = r OFF
out_mode_05	1	Start the circulator
out_sp_01	xxx.x	Set working temperature  *
out_sp_02	xxx.x	Set high temperature warning limit 
out_sp_03	xxx.x	Set low temperature warning limit 
in_sp_01	none	Ask for working temperature 
in_sp_02	none	Ask for high temperature warning limit 
in_sp_03	none	Ask for low temperature warning limit 
in_pv_00	none	Ask for actual bath temperature
in_pv_01	none	Ask for the heater wattage being used

25.4. Status messages

Message	Description
00 MANUAL STOP	Circulator in "OFF" state
01 MANUAL START	Circulator in keypad control mode
02 REMOTE STOP	Circulator in "r OFF" state
04 REMOTE START	Circulator in remote control mode

(* see "Note" on page 18)

25.5. Error messages

Message	Description
-01 TEMP / LEVEL ALARM	Safety temperature or low liquid level alarm
-03 EXCESS TEMPERATURE WARNING	High temperature warning 
-04 LOW TEMPERATURE WARNING	Low temperature warning 
-05 TEMPERATURE MEASUREMENT ALARM	Error in measuring system
-07 I ² C-BUS WRITE ERROR -07 I ² C-BUS READ ERROR -07 I ² C-BUS READ/WRITE ERROR	Internal error
-08 INVALID COMMAND	Invalid command
-10 VALUE TOO SMALL	Entered value too small
-11 VALUE TOO LARGE	Entered value too large
-12 WARNING: VALUE EXCEEDS TEMPERATURE LIMITS	Value lies outside the adjusted range for the high and low temperature warning limits. But value is stored.
-13 COMMAND NOT ALLOWED IN CURRENT OPERATING MODE	Invalid command in current operating mode

26. Cleaning the unit

Before cleaning the unit, disconnect the power plug from the mains socket!

For cleaning the bath tank and the immersed parts of the circulator, use low surface tension water (e.g., soap suds). Clean the outside of the unit using a wet cloth and low surface tension water.

Prevent humidity from entering into the circulator CT 52 E.

27. Maintenance

The transparent-thermostat CT 52 is designed for continuous operation under normal conditions. Periodic maintenance is not required.

The tank should be filled only with a bath liquid recommended by SCHOTT Instruments GmbH. To avoid contamination, it is essential to change the bath liquid from time to time..

Repairs

Before asking for a service technician or returning a SCHOTT Instruments GmbH circulator for repair, please contact an authorized service station.

SCHOTT Instruments GmbH
Hattenbergstrasse 10
D-55122 Mainz / Germany

Tel. +49 61 31 / 66 51 11


Fax +49 61 31 / 66 50 01

E-mail: avs@schottinstruments.com

www.schottinstruments.com

When returning a unit, take care of careful and adequate packing. SCHOTT Instruments GmbH is not responsible for damages that might occur from insufficient packing.

28. Disposal of the unit

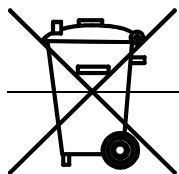Disposal with household waste (unsorted waste) or similar collections of municipal waste is not permitted!

Contact an authorized waste management company in your country.

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